

# Alaska Scientific Crime Detection Laboratory

## Change in Instrument Status Form

Issued: 12/15/2017  
Effective: 12/15/2017

Version: CIISF 2017 R0  
Status: Active

Date: 1-21-2020 Instrument Serial Number: 100346

Instrument Location: Anchorage Jail

Supervisor Name and Agency: Sergeant David Noll (6520) - Anchorage Police Department

### Describe the Status Change or Issue with the Instrument:

On 1-20-2020 Ofc. Kendall (BTO# 9451) was attempting to use Datamaster 100346 at the Anchorage Jail for an OUI breath sample, APD case 20-2394. The touchscreen was not working but Ofc. Kendall was able to tab through options to start a test sequence. The results never printed and Ofc. Kendall was not able to print the last copy. He restarted the Datamaster and the touchscreen still did not work. He tabbed to the Copy Button and printed the last copy which was from a different incident. I responded to the jail and observed the touchscreen was not working. I attempted a hard reset and had the same result. I could not tab to the main menu to look for the report that way and could not initiate a VOC. I selected Copy and observed it printed a prior incident. I also could not use the menu to place the Datamaster Out of Service.

If instrument was placed in service or removed from service, fill out the appropriate sections below.

In Service Date/Time: \_\_\_\_\_

Out of Service Date/Time: 1-21-2020 at 0100 hours

**For Use by SCDL**

**Additional Notes**

---

---

---

---

---

---

---

---

---

---

Email completed form to [dps.scdl.toxicology@alaska.gov](mailto:dps.scdl.toxicology@alaska.gov)

For questions contact the Breath Alcohol Program at 907-269-5740