

Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

Issued: 7/10/2015
Effective: 7/10/2015

Version: DMCIF 2015 R1
Status: Active

Instrument S/N 100417

Received From NPAS

Date 8/3/15

Analyst Colleen O'Bryant

1. Reason for instrument return:

Returned from NPAS

2. Change instrument and case location in JusticeTrax Breath Alcohol Database

3. Ensure change in instrument status form is in JusticeTrax N/A

4. Upload N/A

5. Update agency information and polling status in DM Host Calling Program N/A

6. Check for VOCs N/A

From NPAS

7. NPAS Repair Paperwork (attached)

Work Ord No 49664

National Patent Analytical Systems, Inc.
Service Action Report

Serial No 100417

Date 6/18/15

Customer Alaska Dept of Public

Phone 907-269-5582

Contact Bolz

War Exp 5/4/13

Mx Exp

Ship Damage? Y N

Cust Pick Up?? Y N

Initial Comments: note any shipping damage or estimates required, along with accessories

Describe problem seen and/or attach documents received

blank/white screen, lock up

Keyword

Finding

Error Code

Enter Hours 2.0

Repair Comments and work Performed

Locks up. Replaced SI pcb assy. Touch screen function intermittent/delayed. Replaced Membrane. Inspected unit, fixed tubing, updated software, set voltages, calibrated, ran standard tests ok

Solution Lot No 15040

Voltage Set Values In				Voltage Set Values Out			
Lamp	1.750	Brth Vol	1.530	Lamp	1.970	Brth Vol	1.540
Meter	1.022	Filter Dif	0.383	Meter	-0.355	Filter Dif	0.494
Stability	0.002			Stability	0.002		

Part No	Qty Ord	Pick	Back Ordered	Total shin	Description	Price Ea
37058	1.0				CLIP, NYLON CABLE .625 DIA (BT HOLDER)	0.000
37553	1.0				MEMBRANE,8.4 IN TOUCH REPLACEMENT replacement	0.000
43602	1.0				KIT, SI CONTROLLER PCB W/PIGGYBK & SERIAL replacement	0.000
9DMT	1.0				DATAMASTER DMT SERVICE RETURN 100417	
NOTE	1.0				NANUK CASE	
r500-1	1.0				Service Charge for Repair	0.000
r500-2	1.0				Service Charge for Repair, Additional Hours	0.000

Date Completed 7/9/15

Technician R. MARTIN

Pick up Signature:

Packing List

From: National Patent Analytical Systems Inc.
2090 Harrington Mem. Rd., Mansfield, Ohio 44903

Cust No. AK6056A Ord Date 6/18/15
Sales Order No. 49664

Ship Date 7/14/15
Packing List No. 74286

Cust Purch Ord No. COLLEEN

Ship To:

Alaska Dept of Public Safety
Scientific Crime Detection
4805 Dr. Martin Luther King Dr.

Anchorage Ak 99507

Attn:

Comments Service for 100417

Bill To:

Alaska Dept of Public Safety
Scientific Crime Detection
4805 Dr Martin Luther King Dr.

Anchorage Ak 99507 **Freight**

25.00

Ship Via FedEx

Part / Drawing No / Rev No	Quantity Ordered	This Shipment	Back Ordered	Total Ship To Date	Description	NP Invent Ref
No Draw # / NONE	1.0	1		1.0	CLIP, NYLON CABLE .625 DIA (BT HOLDER)	37058
Serial Nos.						
No Draw # / NONE	1.0	1		1.0	MEMBRANE, 8.4 IN TOUCH REPLACEMENT	37553
replacement						
Serial Nos.						
43602 / A	1.0	1		1.0	KIT, SI CONTROLLER PCB W/PIGGYBK & SERIAL	43602
replacement						
Serial Nos.						
No Draw # / NONE	1.0	1		1.0	DATAMASTER DMT SERVICE RETURN	9DMT
100417						
Serial Nos.						
No Draw # / NONE	1.0	1		1.0	NANUK CASE	NOTE
Serial Nos.						
No Draw # / NONE	1.0	1		1.0	Service Charge for Repair	r500-1
Serial Nos.						
No Draw # / NONE	1.0	1		1.0	Service Charge for Repair, Additional Hours	r500-2
Serial Nos.						