

Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

Issued: 1/13/2017
Effective: 1/17/2017

Version: DMCIF 2017 R0
Status: Active

Instrument S/N 100424

Received From Intoximeters

Date 02/11/21

Analyst Derec Walton

1. Reason for instrument return:

Refurbished by manufacturer for filter wheel errors

2. Change instrument and case location in JusticeTrax Breath Alcohol Database

3. Ensure change in instrument status form is in JusticeTrax

4. Upload

*Instrument was not in the field
DMV 02/11/21*

5. Update agency information and polling status in DM Host Calling Program

6. Check for VOCs

From Intoximeters

7. Intoximeters Repair Paperwork (attached)

SRO Number: SRO-043864

Received Date: 01/14/2021

Completion Date: 02/04/2021

SRO Type: REPAIR

SRO Description: DMT Preventive Maintenance

Customer ID: C000003651

Customer PO: 01142021-Walton

Contact: Derek Walton

Email: derek.walton@alaska.gov

Ship To:

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507

Units on SRO

| | | |
|---|--------|------------|
| 1 | LEGDMT | LEGACY DMT |
|---|--------|------------|

| | | |
|--------------------------|---|---|
| Service Line: 1 | | **All Instruments are Calibrated to Factory Specifications** |
| Unit Type: | (LEGDMT) LEGACY DMT | |
| Serial Number: | 50300315 | |
| Reason: | DMT Status Messages | |
| | FltWl er | |
| Reason Notes: | Preventative Maintenance, filter wheel error | |
| Resolution: | Prev M | |
| | planned | |
| Resolution Notes: | Removed filter wheel assembly for cleaning and lubrication. Cleaned & adjusted filter wheel solenoid. Updated filter wheel solenoid mounting screws. Updated stepper motor adjustment lock down screw. Replaced filter wheel motor pinion. Cleaned, lubricated, & adjusted filter wheel moving parts. Removed, inspected, and applied metallic seal tape to mirror blocks of Sample Chamber. Inspected 5-Way Valve. Performed Calibration/Adjustment, Wet and Dry Gas Accuracy Checks, Breath Tests, and outgoing Diagnostics Test. | |

| Parts: | <u>Qty</u> | <u>Part Number</u> | <u>Part Description</u> |
|--------------------------|--|--------------------|--|
| | 1 | 90-8000-00 | CONTRACT PREV MAINT DMT |
| Reason: | DMT Functional | | |
| | DigPots | | |
| Reason Notes: | Requires Digital Pots Mod. | | |
| Resolution: | Update | | |
| | DMTVR3 | | |
| Resolution Notes: | Removed Control Board and installed TB-30 DMT Digital Pots Mod. All standard factory checks performed. | | |
| | 1 | 001199-00 | DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN |

Order Packing Slip Report

Intoximeters
Experience • Service • Integrity

2/8/2021 1:21:55 PM

Packing Slip: 220935

Pack Date: 2/8/2021

SRO: SRO-043864



From: Intoximeters, Inc
2081 Craig Road
St Louis MO 63146
USA

Customer: C000003651

Ship To: 2

Alaska Department of Public Safety
4805 Dr MLK Jr Ave
Anchorage AK 99507
USA

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507
USA

Customer PO: 01142021-Walton

Ship Via: FDX

FED EX

Weight: 38.00

Delivery Terms : 2D

FOB : ORIGIN

Packages: 1

Customer Ship Account :

| Line Number | Item | UM | Ordered Quantity | Shipment Quantity | Backorder Quantity |
|--|------------|------------|------------------|-------------------|--------------------|
| 1 | LEGDMT | LEGACY DMT | | | 50300315 |
| 1 | 90-8000-00 | EA | | 1.00 | 0.00 |
| CONTRACT PREV MAINT DMT | | | | | |
| 2 | 001199-00 | EA | | 1.00 | 0.00 |
| DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN | | | | | |