

Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

Issued: 1/13/2017
Effective: 1/17/2017

Version: DMCIF 2017 R0
Status: Active

Instrument S/N 100697

Received From Intox

Date 10/11/21

Analyst Brandi Barnett

1. Reason for instrument return:

Return from factory refurbishment

2. Change instrument and case location in JusticeTrax Breath Alcohol Database

3. Ensure change in instrument status form is in JusticeTrax N/A

4. Upload N/A

5. Update agency information and polling status in DM Host Calling Program N/A

6. Check for VOCs N/A

From Intoximeters

7. Intoximeters Repair Paperwork (attached)

SRO Number: SRO-046022

Received Date: 05/17/2021

Completion Date: 09/28/2021

SRO Type: REPAIR

SRO Description: DMT Preventive Maintenance

Customer ID: C000003651

Customer PO: 050521-Walton

Contact: Derek Walton

Email: derek.walton@alaska.gov

Ship To:

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507

Units on SRO

1	LEGDMT	LEGACY DMT
---	--------	------------

Service Line: 1	**All Instruments are Calibrated to Factory Specifications**	
Unit Type:	(LEGDMT) LEGACY DMT	
Serial Number:	50100697	
Reason:	Preventive Maintenance	
	planned	
Reason Notes:	Customer Reported Issue: Loud regulator	
Resolution:	Replaced	
	RPL comp	
Resolution Notes:	Replaced Drygas Regulator Assembly. Removed filter wheel assembly for cleaning and lubrication. Cleaned & adjusted filter wheel solenoid. Updated filter wheel solenoid mounting screws. Updated stepper motor adjustment lock down screw. Replaced filter wheel motor pinion. Cleaned, lubricated, & adjusted filter wheel moving parts. Removed, inspected, and applied metallic seal tape to mirror blocks of Sample Chamber. Replaced discolored tubing and inspected 5-Way Valve. Replaced flow control assembly.	
Parts:	<u>Qty</u>	<u>Part Number</u>
	1	90-8000-00
		CONTRACT PREV MAINT DMT
Reason:	DMT Functional	
	DigPots	
Reason Notes:	Requires Digital Pots Mod.	
Resolution:	Update	
	DMTVR3	
Resolution Notes:	Removed Control Board and installed TB-30 DMT Digital Pots Mod. Optimized detector position. Adjusted Tech Screen settings.	
	1	001217-00
		DMT HRDW GAS MOUNT ASSY W/VALVE
Reason:	DMT Mechanical or Physical	
	Touch Sc	
Reason Notes:	Touch Screen not functioning correctly.	
Resolution:	Replaced	
	RPL comp	
Resolution Notes:	Replaced Touch Screen Panel.	
	1	59-0000-08
		DMT HRDW CABLE TIE HOLDER 30320
	1	001199-00
		DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN
	1	59-0057-00
		DMT HRDW FLOW CONTROL ASSY 41396
	1	59-0101-00
		DMT ELEC MEMBRANE 8.4IN TOUCH REPL 37553

Order Packing Slip Report

Intoximeters
Experience • Service • Integrity

10/5/2021 2:50:25 PM

Packing Slip: 238184

Pack Date: 10/5/2021

SRO: SRO-046022



From: Intoximeters, Inc
2081 Craig Road
St Louis MO 63146
USA

Customer: C000003651

Ship To: 2

Alaska Department of Public Safety
4805 Dr MLK Jr Ave
Anchorage AK 99507
USA

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507
USA

Customer PO: 050521-Walton

Ship Via: FDX

FED EX

Weight: 37.00

Delivery Terms : 2D

Shipping Terms : ORIGIN

Packages: 1

Customer Ship Account :

Line Number	Item	UM	Ordered Quantity	Shipment Quantity	Backorder Quantity
1	LEGDMT	LEGACY DMT			50100697
1	90-8000-00	EA	1.00	1.00	0.00
	CONTRACT PREV MAINT DMT				
2	001217-00	EA	1.00	1.00	0.00
	DMT HRDW GAS MOUNT ASSY W/VALVE				
3	59-0000-08	EA	1.00	1.00	0.00
	DMT HRDW CABLE TIE HOLDER 30320				
4	001199-00	EA	1.00	1.00	0.00
	DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN				
5	59-0057-00	EA	1.00	1.00	0.00
	DMT HRDW FLOW CONTROL ASSY 41396				
6	59-0101-00	EA	1.00	1.00	0.00
	DMT ELEC MEMBRANE 8.4IN TOUCH REPL 37553				

Rec'd 10/07/21
JM